

Preparation of single-cell suspensions from mouse spleen without enzymatic treatment

Contents

1. Description
 - 1.1 Background information
 - 1.2 Reagent and instrument requirements
2. Protocol for the dissociation of mouse spleen

1. Description

1.1 Background information

Single-cell suspensions are a prerequisite for many experiments, for example to achieve the highest possible purity and recovery during cell separations with MACS® Technology. The gentleMACS™ Dissociator provides optimized programs to attain single-cell suspensions from various tissues, for example, mouse spleen. In combination with C Tubes, the gentleMACS Dissociator allows the automated tissue dissociation in a closed system, enabling sterile sample handling. A single sample or two samples can be processed in parallel.

1.2 Reagent and instrument requirements

- gentleMACS Dissociator
- gentleMACS C Tubes (# 130-093-237)
- Pre-Separation Filters (# 130-041-407)
- (Optional) ART 1000 REACH pipet tips (Molecular BioProducts, Inc.) for removal of dissociated material from the closed C Tubes.
- Buffer: Prepare a solution containing phosphate-buffered saline (PBS), pH 7.2, 0.5% bovine serum albumin (BSA), and 2 mM EDTA by diluting MACS BSA Stock Solution (# 130-091-376) 1:20 with autoMACS™ Rinsing Solution (# 130-091-222). Keep buffer cold (2–8 °C).

▲ **Note:** EDTA can be replaced by other supplements such as anticoagulant citrate dextrose formula-A (ACD-A) or citrate phosphate dextrose (CPD). BSA can be replaced by other proteins such as gelatine, mouse serum or fetal bovine serum.

2. Protocol for the dissociation of mouse spleen

▲ For details on the use of the gentleMACS Dissociator, refer to the gentleMACS Dissociator user manual.

▲ For cell culture experiments subsequent to tissue dissociation, all steps should be performed under sterile conditions.

▲ The weight of one mouse spleen amounts to 80–120 mg (female BALB/c mouse, 6–7 weeks old).

1. Transfer mouse spleen into the gentleMACS C Tube containing buffer:
 - 1–2 mouse spleens: 3 mL
 - 3–4 mouse spleens: 6 mL
 - 5–6 mouse spleens: 9 mL

2. Tightly close C Tube and attach it upside down onto the sleeve of the gentleMACS Dissociator.
 - ▲ **Note:** It has to be ensured that the sample material is located in the area of the rotor/stator.
3. Switch on the gentleMACS Dissociator and choose one of the gentleMACS Programs:
 - 1–2 mouse spleens: **m_spleen_01**.
 - 3–6 mouse spleens: **m_spleen_04**.
4. Run the gentleMACS Program **m_spleen_01** or **m_spleen_04**.
5. After termination of the program, detach C Tube from the gentleMACS Dissociator.
6. (Optional) Perform a short centrifugation step to collect the sample material at the bottom of the tube.
7. Resuspend sample and apply the cell suspension to a Pre-Separation Filter placed on a 15 mL tube (1–2 mouse spleens per C Tube) or to an appropriate cell strainer placed on a 50 mL tube (3–6 mouse spleens per C Tube).
 - ▲ **Note:** Dissociated tissue can be removed from the closed C Tube by pipetting through the septum-sealed opening in the center of the cap of the C Tube. Use ART 1000 REACH 1000 µL pipette tips.
8. Wash Pre-Separation Filter with 5 mL of buffer.
9. Discard Pre-Separation Filter and centrifuge cell suspension at 300×g for 10 minutes at room temperature. Aspirate supernatant completely.
10. Resuspend cells with buffer to the required volume for further applications.

All gentleMACS Protocols are available at www.miltenyibiotec.com.

Warranty

The products sold hereunder are warranted only to be free from defects in workmanship and material at the time of delivery to the customer. Miltenyi Biotec GmbH makes no warranty or representation, either expressed or implied, with respect to the fitness of a product for a particular purpose. There are no warranties, expressed or implied, which extend beyond the technical specifications of the products. Miltenyi Biotec GmbH's liability is limited to either replacement of the products or refund of the purchase price. Miltenyi Biotec GmbH is not liable for any property damage, personal injury or economic loss caused by the product.

MACS is a registered trademark and gentleMACS is a trademark of Miltenyi Biotec GmbH.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for diagnostic or therapeutic use.

Copyright © 2010 Miltenyi Biotec GmbH. All rights reserved.